

## **Read User Manual Before Start-up! How to operate I-Care controller!**

1. Make sure the appliance has been in a stable and levelled upright position for at least 1 hour before plugging it in
2. Ensure that the plug is mounted correctly & securely. Check that the voltage connection is in accordance to the rating label inside the cabinet.
3. The installation location should meet the following requirements:
  - Ensure the ambient temperature is within the range of 41.5°F and 95°F.
  - 2.5 inches of space on all sides and above the appliance.
  - No direct sunlight.
  - Allow the appliance to operate at the set point for at least 12 hours before loading.
  - Clean the inside of the cabinet using luke warm water with a mild detergent. Use soft cloth, rinse with clean water and dry thoroughly.
4. Fit distance pieces to the back of the appliance & the anti-slip pieces under the feet.  
NOTE: This does not apply to all models.
5. If equipped with a reference bottle, remember to fill it. I.e. with 50% water and 50% ethanol.
6. Once you turn on the appliance you will be guided through an installation wizard, which gets you set up.
7. It is important to set time & date on the controller during the installation wizard.

**NB! Quick start guide does not replace manual.  
We advise that you always read the manual  
before operating the appliance!**